

## 1.11

### Editing a Batch in Queue

Std. Familiarization Time: 0:05:00


Approximate Video Time: 0:04:00

[Play the Video](#)

MassHunter Basics for Agilent ICP-MS  
Rev.3.0.2 – February 2021  
50

## Queue

- Ribbon > [Home] tab > [Batch] group > [Add to Queue]
- [Skip Warm-up]
- [Plasma Off at End]
- [Vial# at End]



Agilent Technologies

MassHunter Basics for Agilent ICP-MS  
Rev.3.0.2 – February 2021  
51

## Adding or deleting a sample during acquisition

- Select the first empty line in the [Sample List] pane from the Acquisition Queue
- Click [Start Sample List Editing Mode]
- Add/delete samples
- Click [Exit Sample Editing Mode]

## Others

- Reacquiring the previous batch
- Editing an acquired batch and acquire it again