

## 1.2

### Turn On ICP-MS


Std. Familiarization Time: 0:05:00

Approximate Video Time: 0:02:00

[Play the Video](#)

MassHunter Basics for Agilent ICP-MS  
Rev.3.0.2 – February 2021  
11

## Turn On ICP-MS

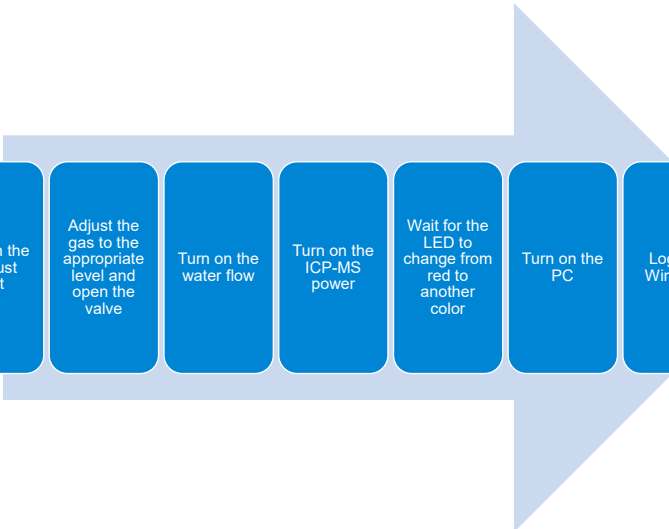

Turn on the exhaust duct

Adjust the gas to the appropriate level and open the valve

Turn on the water flow


Turn on the ICP-MS power

Wait for the LED to change from red to another color

Turn on the PC

Login to Windows

Start MassHunter



Agilent Technologies

MassHunter Basics for Agilent ICP-MS  
Rev.3.0.2 – February 2021  
12