



1.8 Setting up a Batch (3/3) – Sample List

Std. Familiarization Time: 0:05:00
Approximate Video Time: 0:03:30

[Play the Video](#)

MassHunter Basics for Agilent ICP-MS
Rev.3.0.2 – February 2021
36

Sample List

- [Sample Type] (required)
- [Sample Name] (required)
- [Vial#]
- [Level]
- [Use Block List]
- [Validate Method]



Agilent Technologies

MassHunter Basics for Agilent ICP-MS
Rev.3.0.2 – February 2021
37

**Agilent Technologies**

Sample List

Batch: 7900FQ-08_Jan_2021-15_18_403.D
Estimated Time for Batch Acquisition: 1667.870 sec

Skip	Sample Type	Sample Name	Comment	Value	File Name	Replicates	Level	Total DL
1	Sample	Blank		1101				
2	☐ CalBk	Std 1		1102			Level 1	
3	☐ CalStd	Std 2		1103			Level 2	
4	☐ CalStd	Std 3		1104			Level 3	
5	☐ CalStd	Std 4		1105			Level 4	
6	☐ CalStd	Std 5		1106			Level 5	
7	☐ Sample	Blank		1301				
8	☐ Sample	Sample 1		1301				
9	☐ Sample	Sample 2		1302				
10	☐ Sample	Sample 3		1303				
11	☐ Sample	Sample 4		1304				
12	☐ Sample	Sample 5		1305				
13	☐ Sample	Sample 1		1301				
14	☐ Sample	Sample 2		1302				
15	☐ Sample	Sample 3		1303				
16	☐ Sample	Sample 4		1304				
17	☐ Sample	Sample 5		1305				
18	☐							
19	☐							
20	☐							
21	☐							
22	☐							
23	☐							
24	☐							
25	☐							
26	☐							
27	☐							
28	☐							