



1.5

Creating a Batch

Std. Familiarization Time: 0:10:00

Approximate Video Time: 0:06:30

[Play the Video](#)

Batch

The following items are stored in a single batch folder


- Acquisition method (including tuning)
- Data analysis method
- Sample list

Methods for creating a batch

- [Method Wizard]
- [Preset Method]
- [Existing Batch]
- [Blank Template]



Method Wizard

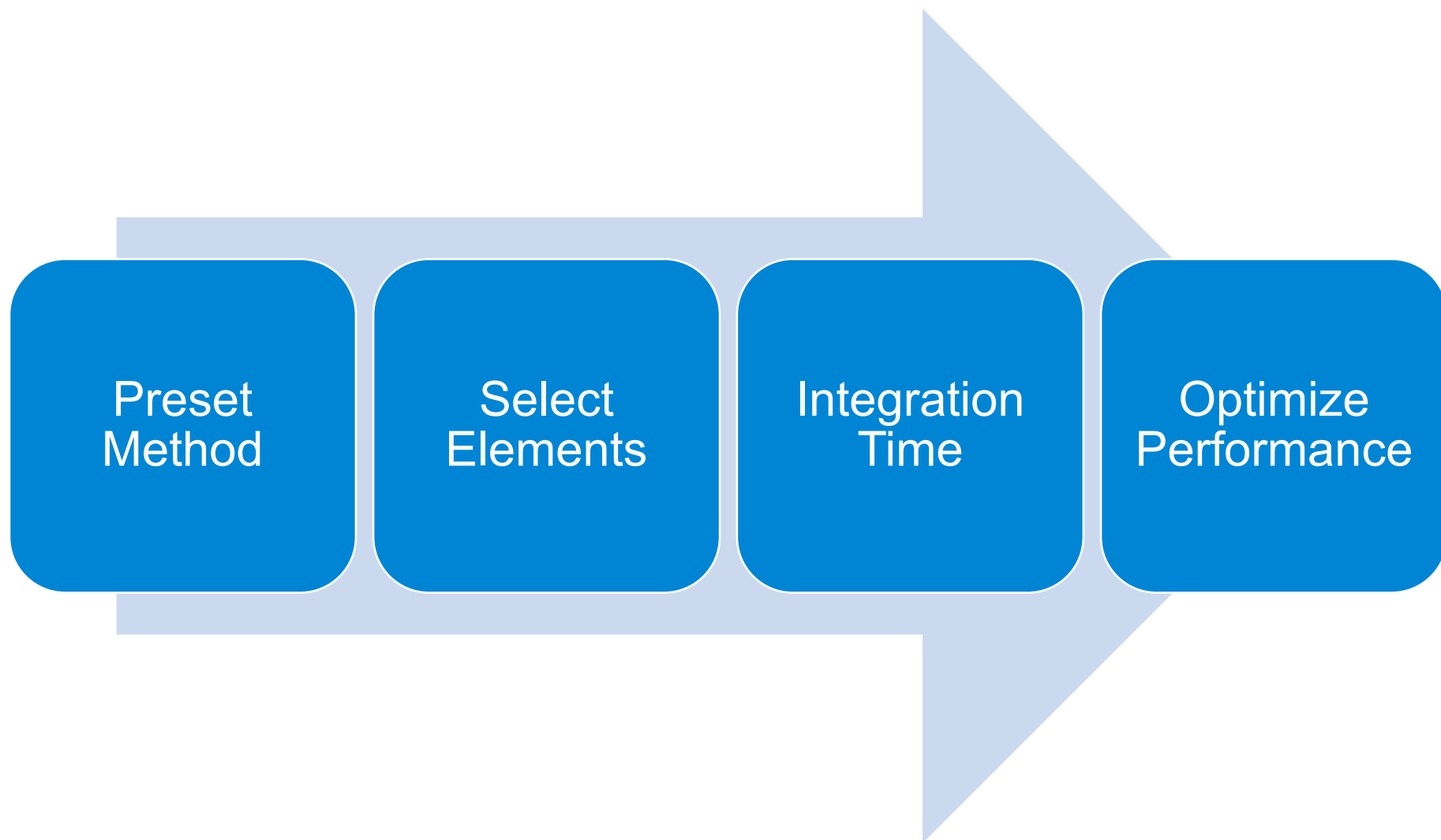

Others

- [File] tab > [Save As Batch Template]
- [Home] tab > [Batch] group > [Create From Template]
- Select the pin of a recent batch template
(This feature has been removed in MassHunter 5.1)
- Support for long file names
- Characters cannot be used for folder or file names