



1.8

Setting up a Batch (3/3) – Sample List

Std. Familiarization Time: 0:05:00

Approximate Video Time: 0:03:30

[Play the Video](#)

Sample List

- [Sample Type] (required)
- [Sample Name] (required)
- [Vial#]
- [Level]

- [Use Block List]

- [Validate Method]

Sample List

[Demo Mode] Online ICP-MS MassHunter - 7900FQ-08_Jan_2021-15_18_40.b

File Home View Report Tools

Nebulizer Pump Speed Autosampler Plasma Configuration Skip Warming up Set as Global Tune New Open Create From Template Save As Add to Queue Advanced Option Resume Pause Stop

Sample Handling Startup Batch Acquisition Queue

Task Navigator

- Hardware
 - Dashboard
- Startup
 - Startup Configuration
 - User Tune Configuration
- Acquisition
 - Setup
 - Tune Modes
 - Element Selection
 - Sample Introduction
 - Monitor
 - Analyte List
 - Calibration
 - SemiQuant
 - QC
 - Worklist Actions
- Sequence
 - Sample List
- Queue
 - Acquisition Queue
- Result
 - Online Data Analysis

Batch - 7900FQ-08_Jan_2021-15_18_40.b

Validate Method Use Block List Import Sample List

Estimated Time for Batch Acquisition: 1667.870 sec

	Skip	Sample Type	Sample Name	Comment	Vial#	File Name	Replicates	Level	Total Dil.
1	<input type="checkbox"/>	Sample	Blank		1101				
2	<input type="checkbox"/>	CalBlk	Std 1		1102			Level 1	
3	<input type="checkbox"/>	CalStd	Std 2		1103			Level 2	
4	<input type="checkbox"/>	CalStd	Std 3		1104			Level 3	
5	<input type="checkbox"/>	CalStd	Std 4		1105			Level 4	
6	<input type="checkbox"/>	CalStd	Std 5		1106			Level 5	
7	<input type="checkbox"/>	Sample	Blank		1201				
8	<input type="checkbox"/>	Sample	Sample 1		1301				
9	<input type="checkbox"/>	Sample	Sample 2		1302				
10	<input type="checkbox"/>	Sample	Sample 3		1303				
11	<input type="checkbox"/>	Sample	Sample 4		1304				
12	<input type="checkbox"/>	Sample	Sample 5		1305				
13	<input type="checkbox"/>	Sample	Sample 1		1301				
14	<input type="checkbox"/>	Sample	Sample 2		1302				
15	<input type="checkbox"/>	Sample	Sample 3		1303				
16	<input type="checkbox"/>	Sample	Sample 4		1304				
17	<input type="checkbox"/>	Sample	Sample 5		1305				
18	<input checked="" type="checkbox"/>								
19	<input checked="" type="checkbox"/>								
20	<input checked="" type="checkbox"/>								
21	<input checked="" type="checkbox"/>								
22	<input checked="" type="checkbox"/>								
23	<input checked="" type="checkbox"/>								
24	<input checked="" type="checkbox"/>								
25	<input checked="" type="checkbox"/>								
26	<input checked="" type="checkbox"/>								
27	<input checked="" type="checkbox"/>								
28	<input checked="" type="checkbox"/>								

Analysis

1/8/2021 4:05:41 PM 3004, Power cycling was detected.

ENG JA